

Planetary Ball Mill ND.4L

Introduction

ND.4L series planetary ball mills could grind or mix solid particles of different granularities and materials, suspensions, and pastes with both dry and wet methods. If a vacuum ball milling tank is used, it will be possible to grind and mix samples in vacuum or inert gases. This lab model provides efficient grinding for superior results up to the nano level, more specifically 30 nm (0.03 μm) according to feedback of users. They are widely used in such fields as geology, metallurgy, soil, building materials, chemical industry, light industry, medicine, electronics, porcelain, battery, environmental protection, and so on.

The planetary Ball Mill performs mixing by continually revolving the sun wheel and rotating the grinding jars concurrently. Both the wheel revolution (centrifugal) speed

and jar rotation (planetary) speed are independently adjustable. This unique mixing capability allows the operator to design precise solutions for various mixing situations. To further improve mixing efficiency, grinding media are added into the grinding jars. Our machines are built in such a way that the strict weight balancing is not required.

Advantages

- Efficient grinding for superior results up to 0.1 μm
- Adjustable speed ratios
- Interval operation function: to prevent the property and quality of the milling materials being affected by overheating, this machine could automatically work in the mode of "operation-shutdown-reoperation".
- Reliable for continuous use
- No strict weight balancing required

TECHNICAL PARAMETERS

Suitable Grinding Jars	50ml, 100ml
Jar Material	Stainless Steel, Stainless Steel-Vacuum Insulated, Agate, Alumina 99.95%, Nylon, Teflon, Tungsten Carbide, Zirconia
Maximum Charge of Ball Milling Tank	Three-fourths of the cubage of the tank (including milling balls)
Feed Size	<5mm; \leq 1mm for crisp materials
Final Fineness	<1 μm ; for colloidal grinding <0.1 μm
Speed Ratio	1:-2 (the grinding jar rotates twice for each sun wheel rotation)
Maximum Rotating Speed	Sun wheel revolution (centrifugal): 300 turns/min \pm 10%; Jar rotation (planetary): 600 turns/min \pm 10%
Operation Model	Mill is controlled by frequency converter and has a total of five operation modes 1) Unilateral operation with shutdown at non-fixed time; 2) Unilateral operation with shutdown at fixed time; 3) Positive and negative alternate operation with shutdown at fixed time; 4) Unilateral interval operation with shutdown at fixed time; 5) Positive and negative alternate interval operation with shutdown at fixed time.
Power	110V, 60HZ, Single-phase, AC, 0.55KW
External Size of Ball Mill	30"x19"x25"/760mmx480mmx640mm
Net Weight	133lb/60KG
Noise Level	\leq 65dB

Torrey Hills Technologies, LLC
6370 Lusk Blvd., Suite F-III
San Diego, CA 92121

Tel (858) 558-6666
Fax (858) 630-3383
Email sales@torreyhillstech.com

